

Megazyme

www.megazyme.com

ASSAY OF
AMYLOGLUCOSIDASE

using

***p*-NITROPHENYL
β-MALTOSE**

plus

Thermostable β-Glucosidase

R-AMGR3 08/18



NOTE:

The AMG assay reagent previously supplied by Megazyme contained *p*-nitrophenyl- β -maltoside (4 mM), plus **almond seed β -glucosidase** (25 U/mL). While this substrate worked well, its use was limited to a narrow pH and temperature range. Also, stability was limited by the purity of the almond β -glucosidase then available. Recently, this substrate has been improved by replacing almond β -glucosidase with a thermostable β -glucosidase. The advantages are:

1. The high purity of the thermostable β -glucosidase means that the substrate is stable for longer periods of time.
2. Because the enzyme is thermostable, the reagent can be used at temperatures up to 60°C (the preferred temperature for the assay of AMG).

However, because thermostable β -glucosidase is unstable at very low salt concentrations, salt and buffer are added to the dry reagent. When dissolved, the substrate solution will have a pH of ~ 6.0 (the optimal pH for stability of the β -glucosidase).

SUBSTRATE:

***p*-Nitrophenyl- β -maltoside (4 mM), plus thermostable β -glucosidase (5 U/mL).**

Dissolve the contents of one vial in 10 mL of distilled water, divide into aliquots of 2-3 mL and store frozen. Store on ice during use.

BUFFER:

100 mM Sodium acetate buffer (pH 4.5).

Add 5.9 mL of glacial acetic acid (1.05 g/mL) to 900 mL of distilled water. Adjust the pH to pH 4.4 by addition of 1 M (4 g/100 mL) NaOH solution. Adjust the volume to 1 L and store in a well sealed bottle at 4°C.

SAMPLE PREPARATION:

Add 1 mL of enzyme preparation to 9 mL of 100 mM sodium acetate buffer (pH 4.5) and mix well. Repeat this dilution step until the enzyme is suitably diluted for assay.

ASSAY PROCEDURE:

1. Pre-equilibrate enzyme solution at 40°C for 5 min.
2. To 0.2 mL of pre-equilibrated substrate solution add 0.2 mL of suitably diluted (and pre-equilibrated) enzyme solution. Mix well and incubate at 40°C for exactly 10 min.

3. Terminate the reaction and develop the colour by adding 3.0 mL of 2% tris base buffer (pH ~ 8.5; Megazyme **B-TRIS500**).
4. Measure the absorbance at 400 nm against a reagent blank.

NOTE:

The reagent blank is prepared by adding 3.0 mL of Tris base buffer (2%) to 0.2 mL of reagent mixture with vigorous stirring, followed by the enzyme solution (0.2 mL) with stirring.

CALCULATION OF ACTIVITY:

$$\text{Activity (U/mL)} = \frac{\Delta A_{400}}{10} \times \frac{3.4}{0.2} \times \frac{1}{18.1} \times \text{Dilution}$$

where:

U = International units of enzyme activity. One Unit is the amount of enzyme which releases one μ mole of *p*-nitrophenol from the substrate per minute at the defined pH and temperature

ΔA_{400} = Absorbance (reaction) - Absorbance (blank)

10 = Incubation time

3.4 = Final reaction volume (mL)

0.2 = Volume of enzyme assayed (mL)

18.1 = E_{mM} *p*-nitrophenol in 2% tris base (pH ~ 8.5) at 400 nm

The Units of amyloglucosidase activity obtained using the above assay, can be related to activity on maltose (10 mg/mL) or soluble starch (10 mg/mL) at 40°C and pH 4.5, using the following equations:

Enzyme Units on Maltose = 1.8 x Units on ***p*NP- β -maltoside**.

Enzyme Units on Starch = 11.5 x Units on ***p*NP- β -maltoside**.

REFERENCE:

McCleary, B. V., Bouhet, F. & Driguez, H. (1991). "Measurement of amyloglucosidase using *p*-nitrophenyl β -maltoside as substrate" *Biotechnology Techniques*, **5**, 255-258.

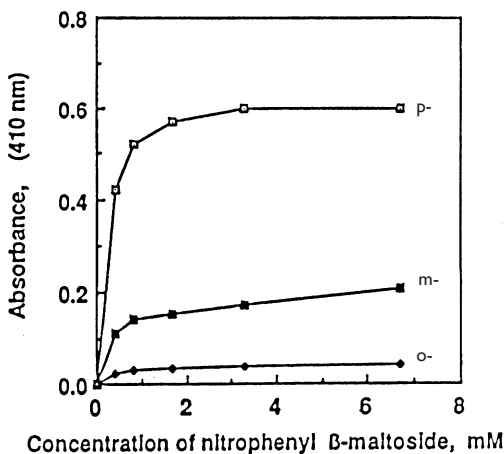


Figure 1. Optimisation of the concentration of nitrophenyl β -maltoside in the reagent mixture.
 Substrates: ■, *m*-; ♦, *o*-; □, *p*-nitrophenyl β -maltosides



**Bray Business Park, Bray,
 Co. Wicklow,
 IRELAND.**

**Telephone: (353.1) 286 1220
 Facsimile: (353.1) 286 1264
 Internet: www.megazyme.com
 E-Mail: info@megazyme.com**

WITHOUT GUARANTEE

The information contained in this booklet is, to the best of our knowledge, true and accurate, but since the conditions of use are beyond our control, no warranty is given or is implied in respect of any recommendation or suggestions which may be made or that any use will not infringe any patents.